

## Supplies Needed for Water Sample Collection



Sterile, clear plastic sample bottle containing Sodium Thiosulfate Powder inside Zip seal bag with absorbent



Refrigerated Cooler with Ice Pack



Outreach Lab Form or Lab Form 507



- Check expiration date on water collection bottle. Samples will not be tested if they do not have an expiration date on the bottle or it is expired.
- <sup>CP</sup> Samples not requiring chain of custody precautions must arrive in the laboratory within 30hrs of collection and are to be kept at <8°C during transport. Use FedEx label provided. Recommend sample collection on Monday, Tuesday, or Wednesday and mail the same day.
- Samples requiring chain of custody precautions must arrive in the laboratory (DLS or another certified laboratory) within 6hrs of collection and are to be kept at <8°C during transport.

Co Pa
East Ball Ball And
Radial Politi Radii Laberatory Ri Barco Bart, Contextuaring Inna, 19 Radia: Data State
Ph Bas 1920 Water Baster Integr
Prost Anno 2000 Water Bucker Indegr
Plan NO 101 Int Aug HO 101 Int
Constant of Advances
manore John Dol
mar by the contraction of a first and the state of the st
Segurituding to
Cary 311/20
Marker The Officer A 2 2
Figst Market of Land Lands Taxabar Same Same Same
Price Indexts of Index actions That a Strating Team Indexts Team Team [1] Index Inc. on a strate [1] under
No. No. 1 Land & the state   Mar.
Dask Ramph Issue Son Rate[] = Parts[] - Cably our sense
(New Actor in a station in the
Cheve Autorities parties are plurible from the first term
nace water row carries
[] (but here if an importantly that all monds have
2 Post Rain Time Drive
A NOT ACAN TANK PUT
Laferatory Radings
1 11
1
1 11
1 1
bether the later being the state of the stat
and a second second second





- Identify sample source, Authorized Collector and Sanitarian ID, Date and Time of Collection, Owner/Occupant, County, and Submitter ID on sample form (LAB 507) to be entered into the Division of Laboratory Services' OUTREACH system at http://chfs.ky.gov/dph/info/lab.
- 2. Select a tap or hose bib which has been in use and does not leak. Do not sample from a drinking fountain or gate valve or at a swivel faucet. Remove all attachments from sample tap prior to sampling.
- 3. Flush tap for 25 minutes before collecting sample. Do not flush tap if source of contamination is suspected to be within the lines of the sampling site.
- 4. Fill the sample bottle to the 100 mL line and close the bottle securely.
- 5. Check the bottle for leaks by shaking and inverting several times.
- 6. Insert water bottle in the zip seal bag with absorbent provided with the water sample kit and seal.
- 7. Place sample bag/s on top of frozen ice pack in cooler and replace styrofoam lid.
- 8. Place printed OUTREACH form on top of styrofoam cooler lid.
- 9. Seal box and place mailing label on top.



Swimming pools and Beaches require 1 (one) 100 mL sample. Testing that will include Pseudomonas, for Public spas, whirlpool, hot tub or other therapeutic pools requires 2 (two)100 mL samples.









If you have questions about submission of water samples, contact the Division of Laboratory Services at (502)-564-4446